


ADPT-OM-ME/W-SG

Adapter between the Omnetic Connectors on the μ PA of the ME-System and on the Headstages of the Wireless-System and the Signal Generator

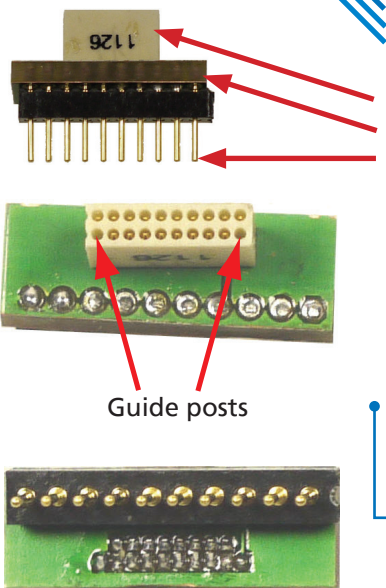

Omnetic socket on W8- or W16 headstage
ADPT-OM-ME/W-SG
Single row precision socket for MEW-SG


Please insert into this Omnetic socket the μ PA or the wireless W8 or W16 headstage, which are equipped with Omnetic connectors.

Important: Take care for the correct orientation!

There are two guide posts on the left and on the right side of the lower pin row of the Omnetic socket, which have to fit to the guide posts of the connector.

Plug the MEW-SG into this single row precision socket.

Important: Take care for the different orientation of the W8 and the W16 headstages!
Please connect the MEW-SG into the ADPT-OM-ME/W-SG as shown on the photo below!

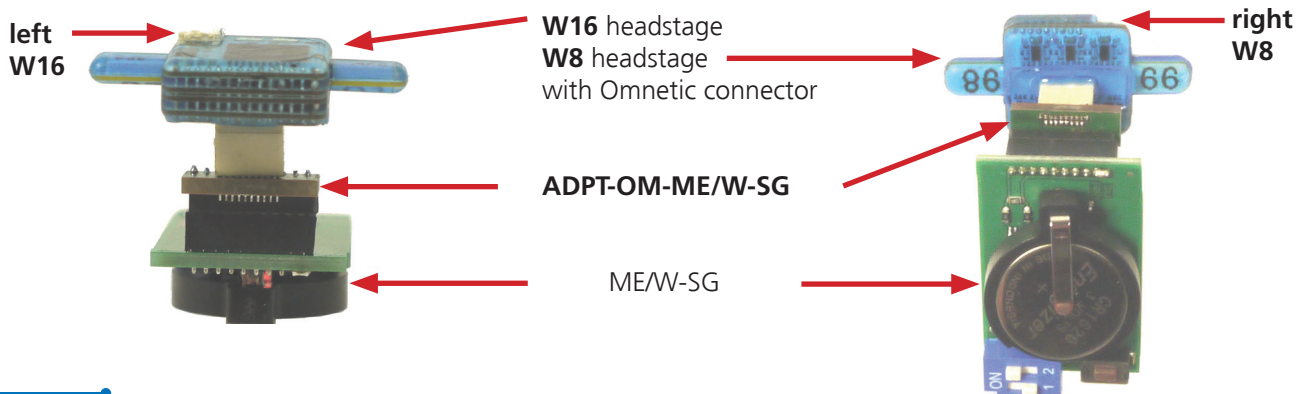

When looking on the service side of the MEW-SG, the white connector for the battery on top of the **W16** headstage has to be on the **left** side!

When looking on the service side of the MEW-SG, the white connector for the battery on top of the **W8** headstage has to be on the **right** side!

Scope of delivery: 1 ADPT-OM-ME/W-SG